

A-0004 Comparator Light Source (CLS)

(For use with C-1001, C-7518, C-7540 and C-7599 only)

1. Prepare the test ampoule according to the kit instructions, then gently place it, flat end first, into the center tube of the comparator.
2. Attach the CLS to the top of the comparator by sliding it in place until it comes to a stop.

Note: A tab-less comparator lid must be in place when the CLS is in use. A spare tab-less lid is enclosed in this package.



3. Press the button on the end of the CLS to turn it on.
4. Hold the comparator assembly up and view from the bottom as illustrated in the kit instructions. To obtain a result, rotate the comparator until the best color match is found.
5. Turn the CLS off and remove it from the comparator.



Simplicity in Water Analysis

www.chemetrics.com
4295 Catlett Road, Midland, VA 22728 U.S.A.
Phone: (800) 356-3072; Fax: (540) 788-4856
E-Mail: orders@chemetrics.com
Oct. 2018, Rev. 3